

Data Sheet

Gravity flow Strep-Tactin® Sepharose® Column

2-1202-001, 2-1202-051, 2-1202-101, Cat. No.:

Lot No.: 1202-

2-1202-550 Version: 11.2 Revision Date: 25.02.2020

IBA Headquarters IBA GmbH

Rudolf-Wissell-Str. 28 37079 Goettingen Germany

Tel. +49 (0) 551-5 06 72-0 Fax +49 (0) 551-5 06 72-181 **IBA US contact information** Fax 1-888-531-6813

E-mail: info@iba-lifesciences.com http://www.iba-lifesciences.com

Description	Ready-to-use column with immobilized streptavidin variant called Strep-Tactin® (5 mg/ml resin) which has been especially optimized for the purification of Strep-tag®II fusion proteins*.
Support	Sepharose 4 FF, 4% agarose
Form	Pre-packed in buffer, pH 8.0: 100 mM Tris-HCl pH 8.0, 1 mM EDTA, 150 mM NaCl, 0.02% sodium azide.
Biotin binding activity	> 300 nmol/ml resin
Stability	6 months after shipping
Storage	recommended: 2- 8 °C
Shipping	room temperature
Hazards	Product is not classified as hazardous according to (EC) No 1272/2008 [CLP]. A Material Safety Data Sheet is provided.

Elution	Strep-Tactin® Elution Buffer with Desthiobiotin (Buffer E), pH 8.0: 100 mM Tris-HCl, 150 mM NaCl, 1 mM EDTA, 2.5 mM desthiobiotin It may be advantageous to use 5-10 mM desthiobiotin to get the target protein eluted at higher concentration.
Regeneration	Strep-Tactin® Regeneration Buffer with HABA; Buffer R If HABA cannot be efficiently removed from Strep-Tactin Sepharose by using Buffer W, we recommend using Buffer W at pH 10.5 (or alternatively 100 mM Tris base) for efficient removal of HABA.

Voss, S. & Skerra, A. (1997) Mutagenesis of a flexible loop in streptavidin leads to higher affinity for the Strep-tag II peptide and improved performance in recombinant protein purification. Protein Eng. 10, 975-982.



Go digital and help the environment. Please download all up-to-date manuals, protocols and other material from http://www.iba-lifesciences.com.

For research use only

Important licensing information

This product is based on Strep-Tactin technology covered by intellectual property (IP) rights and on completion of the sale IBA grants respective Limited Use Label Licenses to purchaser. IP rights and Limited Use Label Licenses for said technology are further described and identified at http://www.ibalifesciences.com/patents.htmlor upon inquiry at info@iba-lifesciences.com or at IBA GmbH, Rudolf-Wissell-Str. 28, 37079 Goettingen, Germany. By use of this product the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses.

Trademark information