

Data Sheet

Gravity flow Protein A Agarose Column – 0.5 mL

Cat. No.: 6-2015-001

Lot No.: 2015-

Version: 2.0
Revision Date: 07.01.2019

Description	Ready-to-use column with Protein A Agarose resin for the purification of monoclonal- and polyclonal antibodies.
Column volume	0.5 mL
Target substance	Polyclonal and monoclonal antibodies (IgG), bound via the Fc-region
Form	Pre-packed in buffer, pH 7.4: PBS pH 7.4, 1 mM EDTA, 0.01 % sodium azide.
Matrix	Rigid, highly cross-linked agarose
Agarose Average particle size ¹(Dv50)	45 µm
Ligand	Alkaline stable recombinant protein A expressed in <i>E. coli</i> using animal-free medium
Dynamic binding capacity ²(DBC10%)	≥70 mg hu IgG/mL resin at residence time 6 min ≥50 mg hu IgG/mL resin at residence time 2.4 min ≥25 mg hu IgG/mL resin at residence time 1 min
Recommended pH working range	2 to 13 (normal working conditions)
Cleaning-In Place stability	No significant change in performance after 100 cycles with 0,5M NaOH (1cycle= 15 min contact time) use at room temperature
Chemical stability	Stable to commonly used aqueous buffers in Protein A chromatography
Storage	recommended: 2- 8°C
Shipping	room temperature
Hazards	Product is not classified as hazardous according to (EC) No 1272/2008 [CLP]. A Material Safety Data Sheet is provided.

¹Median particle size of the cumulative volume distribution.

²Dynamic binding capacity was determined with 1 mg/ml monoclonal hu IgG1 using a 1 ml column at 1, 2.4 and 6 minutes residence time.



Go digital and help the environment. Please download all up-to-date manuals, protocols and other material from <https://www.iba-lifesciences.com/download-area.html>.

For research use only

Trademark information

The owners of trademarks marked by “®” or “TM” are identified at <http://www.iba-lifesciences.com/patents.html>. Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

